


Quick Reference Guide for Using BD FACSuite™ Clinical and BD FACSuite™ Software with the BD FACSLink™ LIS Interface Solution

This guide contains instructions for sending results to a Laboratory Information System (LIS) using the BD FACSLink™ interface embedded in BD FACSuite™ and BD FACSuite™ Clinical software.

Workflow Overview

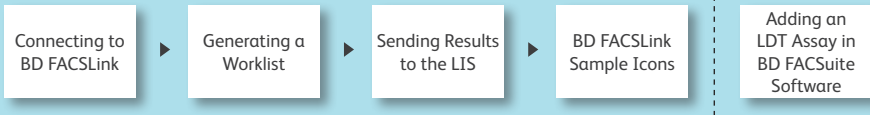

Connecting to BD FACSLink

- 1 Open the software and log in as an administrator.
- 2 Select **Tools > Preferences**, then select the **Worklist** tab.
- 3 Enter the following information:


BD FACSLink Connection

Username: IM_ADMIN

Password: []

TCP/IP Address: http:\\10.10.115.67 Port: 10006 [Test Connection]

- 4 Click **Test Connection** after the information has been entered.
 - A dialog opens indicating that the connection was successful.
- 5 Click **OK** and close the **Preferences** dialog.

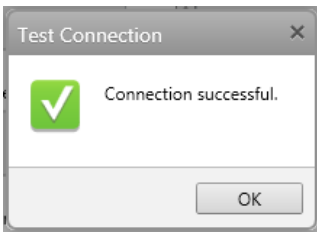

- 6 You will see that the software is connected to BD FACSLink on the status bar and under System Status.




BD FACSuite Clinical Software is CE marked. For In Vitro Diagnostic Use.
BD FACSuite Software is for Research Use Only. Not for use in diagnostic or therapeutic procedures.

© 2017 BD. BD, the BD Logo and all other trademarks are property of Becton, Dickinson and Company.

23-19445-00

Generating a Worklist


- 1 Navigate to the **Worklist** workspace, and create a new worklist.
- 2 Scan barcodes or manually type in the Sample ID column and press Enter.
 - The software will automatically query BD FACSLink for any test orders associated with the Sample ID.
 - Associated patient information from the LIS is added to the worklist entries.
- 3 Continue until the worklist is complete.

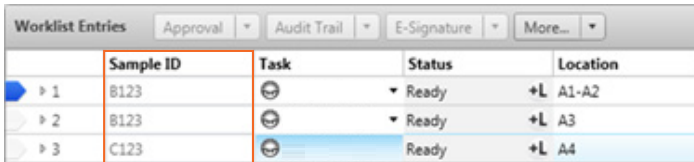

The screenshot shows a table titled "Worklist Entries" with columns: Sample ID, Task, Status, and Location. There are three rows of data. The first row has Sample ID B123, Task (with a smiley face icon), Status Ready, and Location A1-A2. The second row has Sample ID B123, Task (with a smiley face icon), Status Ready, and Location A3. The third row has Sample ID C123, Task (with a smiley face icon), Status Ready, and Location A4. The table has a grey header and a light blue background for the data rows.

Sample ID	Task	Status	Location
B123	☺	Ready	A1-A2
B123	☺	Ready	A3
C123	☺	Ready	A4

- 4 If multiple tests are ordered, a dialog will open to let you select the test to add. Use the **Add All** or **Add Selected** buttons, or double-click one test.


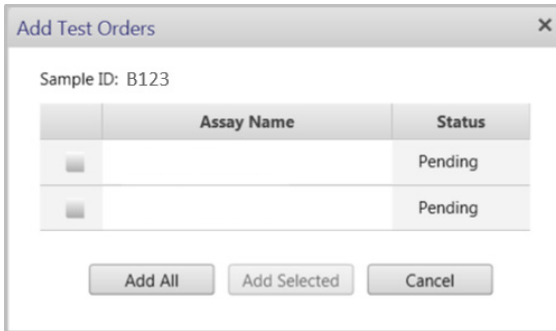


The screenshot shows a dialog box titled "Add Test Orders" with a close button (X) in the top right corner. The dialog contains the text "Sample ID: B123" and a table with two columns: "Assay Name" and "Status". There are two rows in the table, both with "Pending" in the Status column. Below the table are three buttons: "Add All", "Add Selected", and "Cancel".

Assay Name	Status
	Pending
	Pending

Sending Results to the LIS

- 1 After acquisition, analyze the entry and mark the entry **Approved** in either the **Entry Details** panel or the **Worklist Entries** panel. The results are automatically sent to the LIS upon approval of an entry.

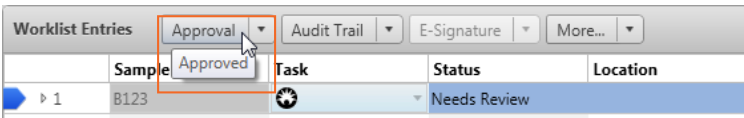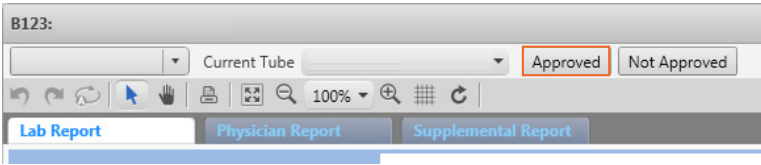

- 2 Select **Tools > BD FACSLink Test Results History** to view information about when the results were transferred for all worklists.


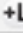
A screenshot of a window titled 'BD FACSLink Test Results History'. It contains a table with the following data:

Status	Worklist Name	Sample ID	Task Name	Approved Time	Delivered Time
Sent	Worklist_001	B123		10/21/2016 11:51:2	10/21/2016 11:51:23 AM
Sent	Worklist_001	C123		10/21/2016 11:51:2	10/21/2016 11:51:22 AM
Sent	Worklist_001	B123		10/21/2016 11:51:1	10/21/2016 11:51:20 AM
Sent	Worklist_001	B123		10/21/2016 11:46:3	10/21/2016 11:46:31 AM
Sent	Worklist_002	123		09/13/2016 12:29:5	09/13/2016 12:29:54 PM



At the bottom right of the window, there are three buttons: 'Refresh', 'Delete', and 'Close'.

BD FACSLink Sample Icons



Icons indicate that an entry was created from a test order.

Worklist Entries			
Approval ▼			
Audit Trail ▼			
E-Signature ▼			
More			
	Sample ID	Task	Status
▶ 1	B123		Ready 

Results are released upon approval of an entry, and icons indicate when results have successfully been sent to BD FACSLink.

Worklist Entries			
Approval ▼			
Audit Trail ▼			
E-Signature ▼			
More			
	Sample ID	Task	Status
▶ 1	B123		Approved 

Icons indicate if results are not sent to BD FACSLink due to a connection problem.

Worklist Entries			
Approval ▼			
Audit Trail ▼			
E-Signature ▼			
More			
	Sample ID	Task	Status
▶ 1	B123		Approved 

Adding an LDT Assay in BD FACSuite Software


To use BD FACSLiNK in BD FACSuite software, you must create an assay with the following characteristics:

- Contains a Statistics View with the Include In Auto Export property turned on.
- Contains LIS Assay Names and Results Names.
- Is a shared assay.


- 1 Select the **Library** icon to go to your library.



- 2 Expand **Assays**, select **User-Defined**, then select your assay.

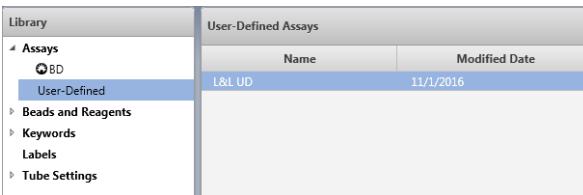


- 3 Select the **Send to LIS** tab and click **Edit**.

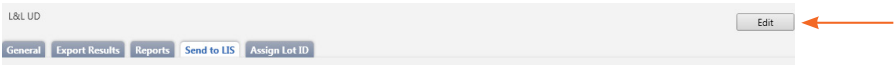

- 4 Enter a BD FACSLiNK Assay Name.

- Only letters, numbers, and underscores are allowed.

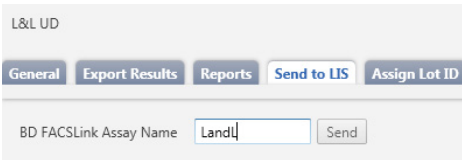

- 5 Check the statistics that you would like imported to your LIS.

- 6 Enter the required names in the **BD FACSLiNK Result Name** column.



A confirmation dialog will be displayed.

BD FACSuite Clinical Software is CE marked. For In Vitro Diagnostic Use.
BD FACSuite Software is for Research Use Only. Not for use in diagnostic or therapeutic procedures.

© 2017 BD. BD, the BD Logo and all other trademarks are property of Becton, Dickinson and Company.
23-19445-00



BD Life Sciences
Biosciences
bdbiosciences.com